MIRO ALTITUDE


Tøuchscreen, single-channel, laser power & energy meter

New Product





CONNECTIVITY

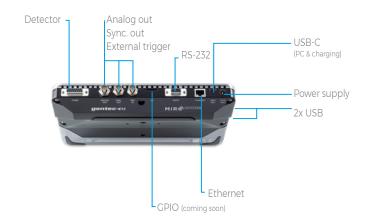

KEY FEATURES

> READS ALL HEADS

Power: thermopiles, photodetectors and pyroelectrics Energy: thermopiles (in SSE mode) and pyroelectrics

LARGE TOUCHSCREEN DISPLAY

10in diagonal 1280 x 800 resolution Touchscreen controls

> INTUITIVE USER INTERFACE

Easy to navigate interface, with 3 display modes: scope, needle and bar chart.
Instant access to the detector settings

REAL-TIME STATISTICAL FUNCTIONS

Max, min, average, standard deviation, RMS and PTP stability, and repetition rate

> MULTIPLE OUTPUTS

Multiple USB ports for computer connection and charging (Ix USB-C, 2x USB-A), BNC analog output, RS-232, Ethernet, programmable I/O (coming soon)

ACCESSORIES



Additional 12V power supply



Power cord extension



USB-C wall charger (US only)



Extra USB-C to USB-A



Pelican carrying case



STAND-R-443



Extra carrying sleeve

145


MIRO ALTITUDE Specifications





	MIRO ALTITUDE
DETECTOR TYPES	Thermopiles, pyroelectrics, photodetectors
DISPLAY	10" high-resolution, anti-glare, touchscreen
POWER METER SPECIFICATIONS	
Power range	4 pW to 150 kW
Meter accuracy	$\pm 0.5\% \pm 3~\mu V$ from 20% to full scale
Statistics	Current value, max, min, average, standard deviation, RMS & PTP stability, time
ENERGY METER SPECIFICATIONS	
Energy range	2 fJ to 30 kJ
Meter accuracy	$1.0\% \pm 50~\mu V$ (< $500~Hz$) $2.0\% \pm 50~\mu V$ (500 Hz to 10 kHz)
Software trigger level	0.1 to 99.9%, 0.1% resolution, default 2%
Repetition rate	10 kHz for data acquisition in real time with time stamp, no missing point
Statistics	Current value, max, min, average, std dev., RMS & PTP stability, pulse #, rep. rate and average power
DETECTOR COMPATIBILITY	
Thermopile	Average power & single shot energy (UP, XLP & HP series)
Photodetector	Average power (PH series)
Pyroelectric	Average power & pulse energy (QE series, except QE8)
GENERAL SPECIFICATIONS	
Digital display size	10.1-inch diagonal LCD - 1280 x 800 pixels
Outputs	Analog out, 0 - 5 V (BNC) Sync out (BNC) RS-232 (DB9) Ethernet (RJ45) USB-C 2x USB-A
Rising edge external trigger	3.3-24 V (BNC)
Serial commands via	USB-C, RS-232 or Ethernet
Data storage via	Internal memory or USB key
Battery type	Rechargeable Li-ion cell
Battery life	6 hours
External power supply	12 VDC power supply included, or UBS-C (min 18 W)
PHYSICAL CHARACTERISTICS	
Mounting holes	1/4"-20 and 2x10-32 threaded holes
Dimensions	268W x 196H x 36D mm
Weight	1.36 kg
ORDERING INFORMATION	
Compatible stand	STAND-R-443
Product page	

MIRO ALTITUDE Main screen

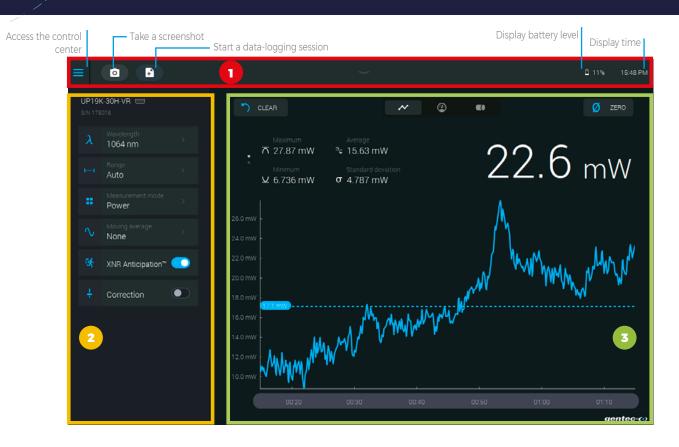

1 NAVIGATION BAR

The upper part of the screen includes a direct access to the control center, data acquisition buttons and various indicators (battery level and time).

2 MEASUREMENT SETTINGS PANEL

Use the various measurement settings available for your detector to set everything related to your measurement.



Wavelength: Enter your wavelength or choose from a list of recently used wavelengths



Range: Set the measuring range to autoscale or to one of the standard values



Measurement mode: Choose what you want to measure: power, SSE, moving average, etc.



Moving average: Choose the desired moving average to use to plot the chart



Trigger: Enter the desired trigger level or choose from a list of recently used values



XNR Anticipation™: Toggle on to measure up to 10x-20x faster without losing any significant accuracy in your readings



Correction: Set a multiplier and an offset value for your measurements



Attenuator: Toggle when using a Gentec-EO calibrated attenuator with your detector

3 DISPLAY AREA

The top part of the display area is the same for all three display modes.



Clear: Use this button to reset the statistics and erase the scope graph's









Zero: Set the current measured value to zero


147

MIRO ALTITUDE Display modes

SCOPE DISPLAY

With this display mode, you can travel in time using the time line at the bottom to view measurements at any point in time while MIRO ALTITUDE continues to measure.


The dotted blue line shows the average value.

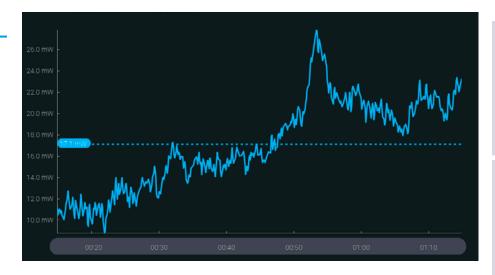

NEEDLE DISPLAY

Faster than an analog needle thanks to XNR Anticipation™! This mode is particularly useful when tuning a laser. The real-time value and statistics are always visible at the top of the screen.


Arrows indicate the minimum and maximum measured values since the last reset. The zoom function sets these values as full scale of the digital gauge.

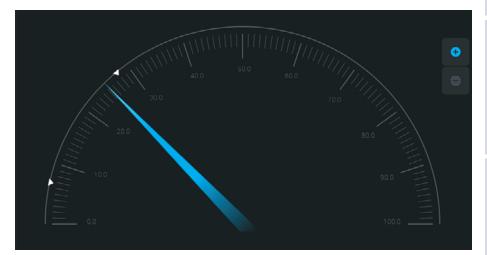

BAR DISPLAY

This is the simplest display mode. Its main advantage is that the current measured value is displayed in huge size, allowing you to read the measurement from a good distance.

Arrows indicate the minimum and maximum measured values.

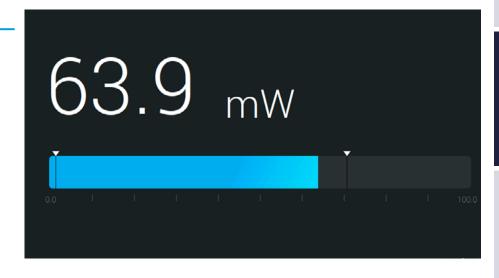

RO ALTITUDE Settings & controls


BUILT-IN FILE MANAGER AND DATA VIEWER

MIRO's built-in file manager lets you access and organize all your screenshots and recorded measurement sessions. You can also copy files on your USB key.



Visualize a recorded measurement session with our built-in data viewer. Data will be displayed in the scope chart display.

There is also a built-in image viewer so you can view your screenshots directly on your MIRO ALTITUDE.

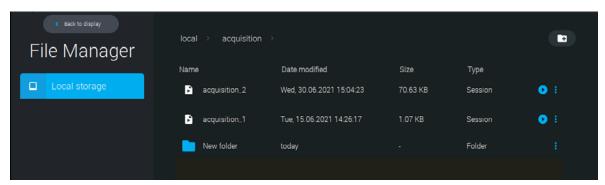

CONTROL CENTER

The control center is accessible from all screens in the top left corner.

Easily navigate between the main screens of the app:



Display




File manager



Settings for the device.

Connect/disconnect your Gentec-EO detector to MIRO ALTITUDE

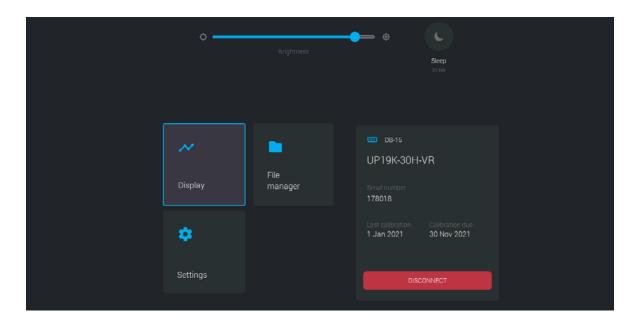


149

MIRO ALTITUDE



DEVICE SETTINGS

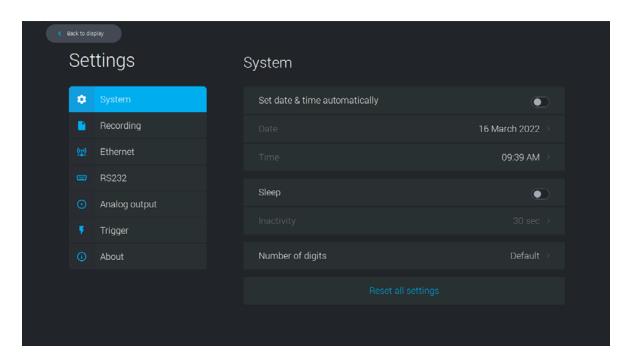



System: Set device settings: language, date, time, sleep, number of digits, etc.



Recording: Set your default recording parameters for power/energy measurement and destination.



Ethernet: Set your Ethernet parameters or let MIRO manage this automatically.



RS-232: Set your RS-232 parameters.



Analog output: Set your analog output parameters. (coming soon)



Trigger: Use an external trigger and set your trigger level.



About: View important information about your device (serial number, firmware version, software version, calibration date) and find support.